



Installing an Eyepiece Reticle

Step 1:
Remove the eyepiece from the microscope's binocular eyepiece tube.



Step 2:
Separate the eyepiece into two pieces by carefully turning the ring at the base of the eyepiece counter clockwise.

Step 3:
Hold the ring with the ledge side up and carefully place the reticle on the ledge. Keep in mind that Nikon reticles should be mounted with the scale side of the glass facing the ledge. Reassemble the eyepiece.



Step 4
Place the eyepiece back into the microscope's binocular tube.

ret·i·cle

noun
(rĕt'ī-kĕl)

An optical element located at the focal plane of an imaging system, containing marks that assist in making measurements of the object under study or help in pointing an instrument. The marks may be as simple as a pair of crossed lines or they may be a complex pattern.*

